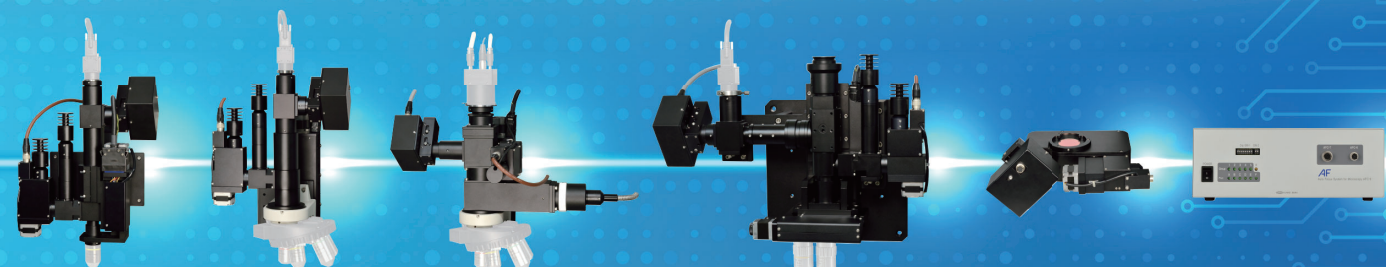
Auto Focus

AUTO FOCUS MICROSCOPE SYSTEM

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Demo room

Visit our demo room to see our demonstration in Tokyo and Osaka office.
Please contact us for more information.

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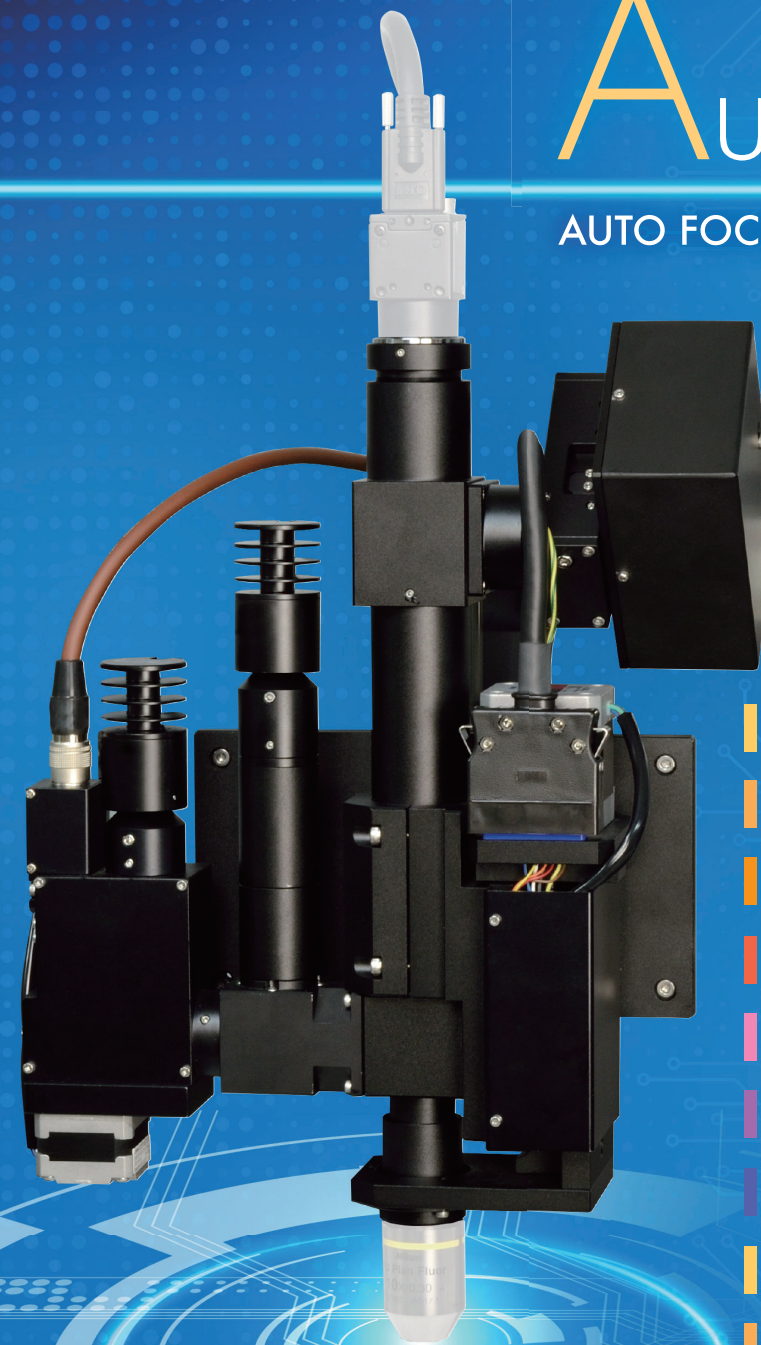
2-line sensor autofocus system achieves superior tracking performance.

Shorter cycle time. Higher productivity.

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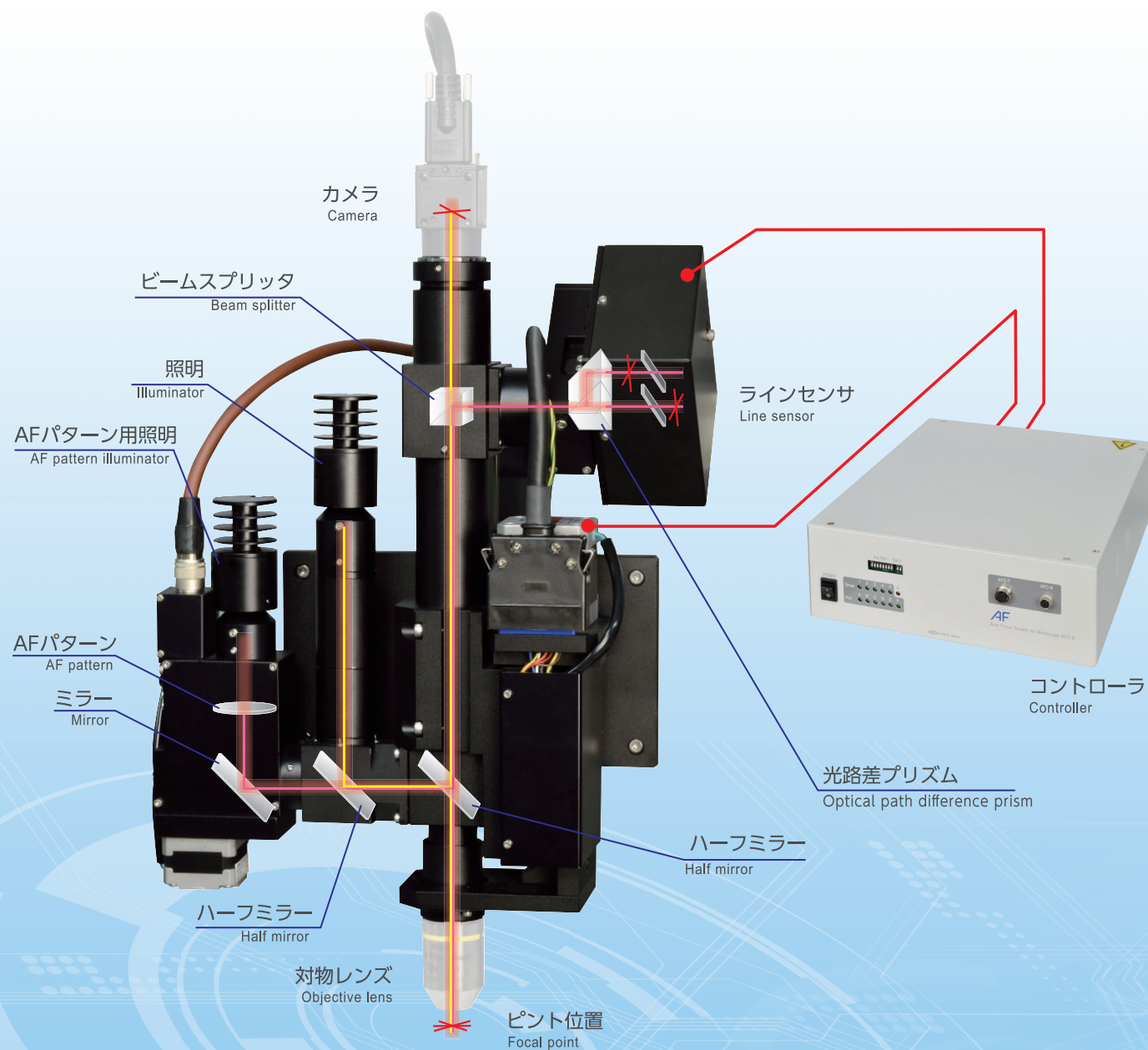
About autofocus microscopes (AF microscopes)

When people hear "autofocus", many may think of the autofocus in digital cameras. Even if a photographer or a subject moves slightly, a digital camera can focus on an image. This does not mean that the camera completely tracks the subject with the autofocus. It is just that the focus area is large, so the image can be focused. However, this is not the case with a microscope. If an objective lens of a microscope is 100x (the total magnification exceeds 1000x), the focus depth (range where the subject comes into focus) is approx. $0.3\mu\text{m}$ (0.0003mm).

This means that if the microscope and the sample are misaligned by just $0.3\mu\text{m}$, the subject is no longer in focus. Our autofocus microscope is a high-technology product that allows for automatic ultramicroscopic focusing adjustment and tracking to keep a moving sample in focus.

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World's leading tracking performance.

World's leading tracking performance. 2-line-sensor method for high-speed tracking of undulations and slopes can maintain continuous focus.

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World's only line-sensor autofocus system.

The world's only autofocus microscope system that utilizes line sensor (one-dimensional image sensors).

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World's leading focusing performance than ever.

Supports both high-reflectance metal surfaces and low-reflectance surfaces such as paper or filters. Also performs well on glass and transparent plastic, which is not possible with other methods. Our autofocus microscope system is high-functioning, and there are few unsuitable samples.

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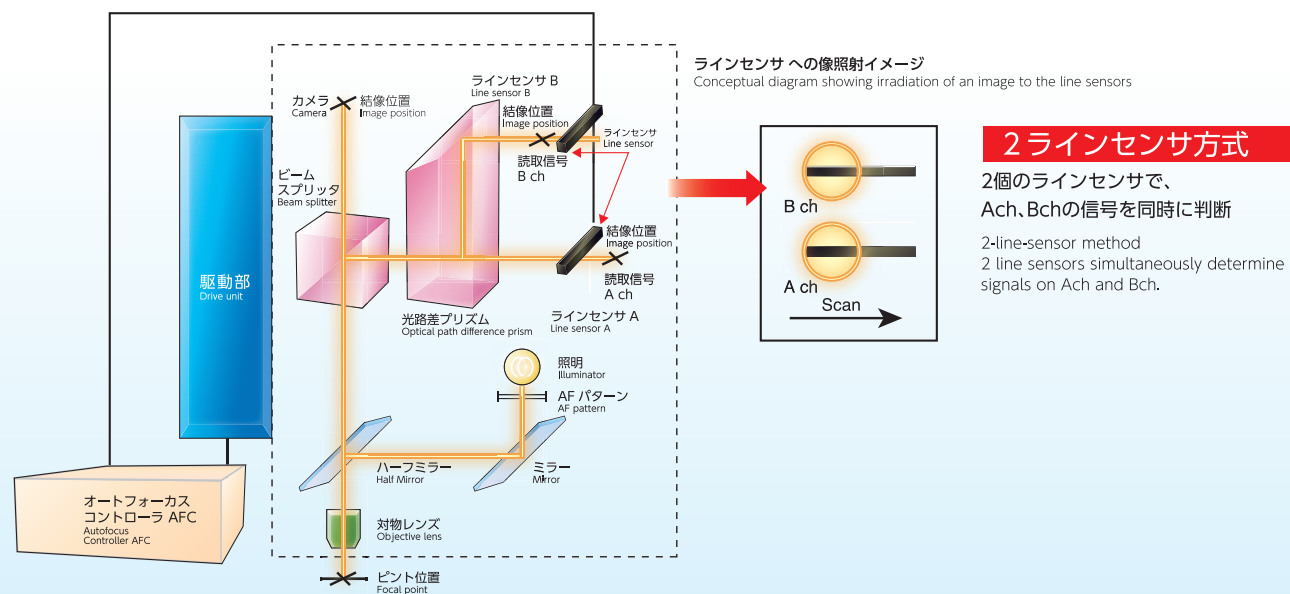
From macro to micro.

Supports a broad range of magnifications from macro to micro.

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Simple idea, huge benefits! 2-line sensor method

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■ Description of the autofocus system

Our autofocus microscope system is mainly composed of a line sensor unit, AF pattern unit, drive unit, controller unit, and lens barrel unit. An AF pattern, which is used for autofocus, is projected on the sample by using the AF pattern irradiation optical system. The AF pattern projected on the sample is split into two optical paths at the optical path difference prism. Light from two images (AF patterns) are received with line sensors. The image signal that is far from the optical path difference prism is called the Ach signal, and the image signal that is near the prism is called the Bch signal. The detection system, including the line sensors, is designed to focus where the Ach and Bch signals are the same. The AF controller performs high-speed signal processing, such as contrast detection, for the Ach and Bch signals obtained from the line sensors, so that the drive unit optimally moves to focus where both signals approximate to focus.

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- ① AF-IZ is designed to move only objective lens unit for autofocus with unified drive unit and microscope tube, to deliver the space saving.
- ② A compact microscope that can be built into an online machining/measuring apparatus.
- ③ This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
- ④ Enables to autofocus on transparent surfaces such as glass and plastics.
- ⑤ Applicable also to surfaces containing potholes or holes.
- ⑥ Supports a broad range of magnifications from macro to micro.
- ⑦ The sensor unit (1-sensor or 2-sensor), AF pattern unit (travel device or without travel device) and illumination unit that can be chosen according to application.
- ⑧ The AF light source (visible or infrared) and the AF pattern unit that can be chosen depending on observation method.

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Objective lenses	Nikon, Olympus, Mitutoyo
Camera	C mount camera (Device size 2/3 inch or less)
Drive unit	Model AF-361ZA
	Limiting resolution 0.078 μm
	Travel Standard 4 mm (changeable up to 10 mm)
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)
Focusing	Line sensor method
Focusing conditions (reference value)	Search range ±250 μm (objective lens 20×)
	Search time 0.3 sec
AF Accuracy	Depth of Focus x 1/4

※ Maximum cable length 15 m.
 ※ Inquire for custom specifications.
 ※ For custom specifications, provide us with your objective lenses and camera.
 ※ The autofocus accuracy and search time vary depending on the sample conditions, etc.

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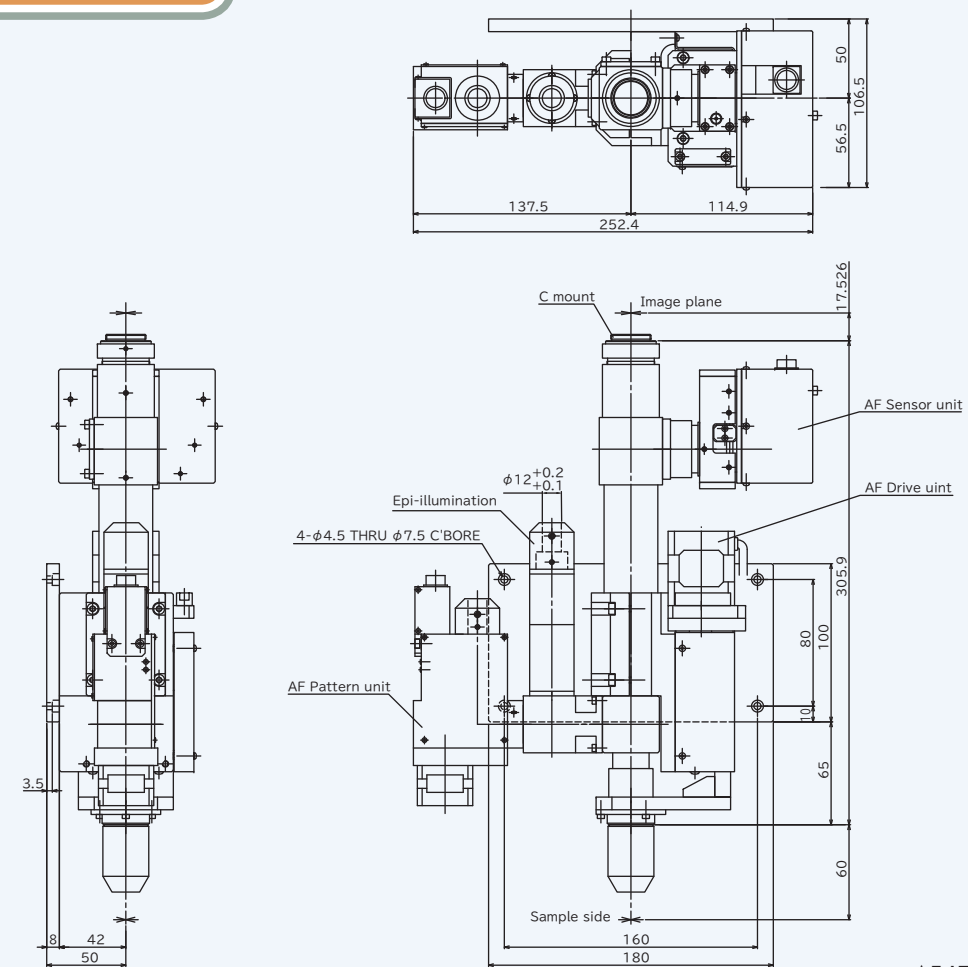
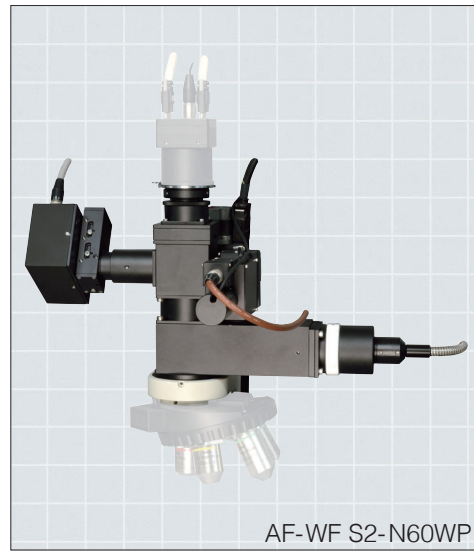
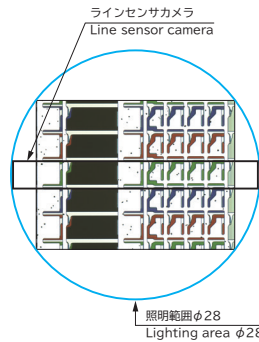
- ① AF-WF is designed to comply with up to φ28 mm or 1.1 inches sensor size, such as C mount area camera and F mount area camera.
 - ② This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
 - ③ Enables to autofocus on transparent surfaces such as glass and plastics.
 - ④ Applicable also to surfaces containing potholes or holes.
 - ⑤ Supports a broad range of magnifications from macro to micro.
 - ⑥ C mount camera or F mount camera can be used with the dedicated camera mount.
 - ⑦ Observation methods (Brightfield, Polarizing and Differential Interference Contrast) that can be chosen depending on the sample.
- *The imaging area vary depending on objective lens and the sample conditions.

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Objective lenses	Nikon, Olympus, Mitutoyo
Camera	C mount camera (Device size 1 inch or less), F mount camera (φ28 mm, equivalent to 1.1 inches)
Drive unit	Model AF-61ZA
	Limiting resolution 0.078 μm
Travel	Standard 4 mm [changeable up to 10 mm]
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)
Focusing	Line sensor method
Focusing conditions (reference value)	Search range ±250 μm [objective lens 20×]
	Search time 0.3 sec
AF Accuracy	Depth of Focus x 1/4

※Maximum cable length 15 m.
 ※Inquire for custom specifications.
 ※For custom specifications, provide us with your objective lenses and camera.
 ※The autofocus accuracy and search time vary depending on the sample conditions, etc.

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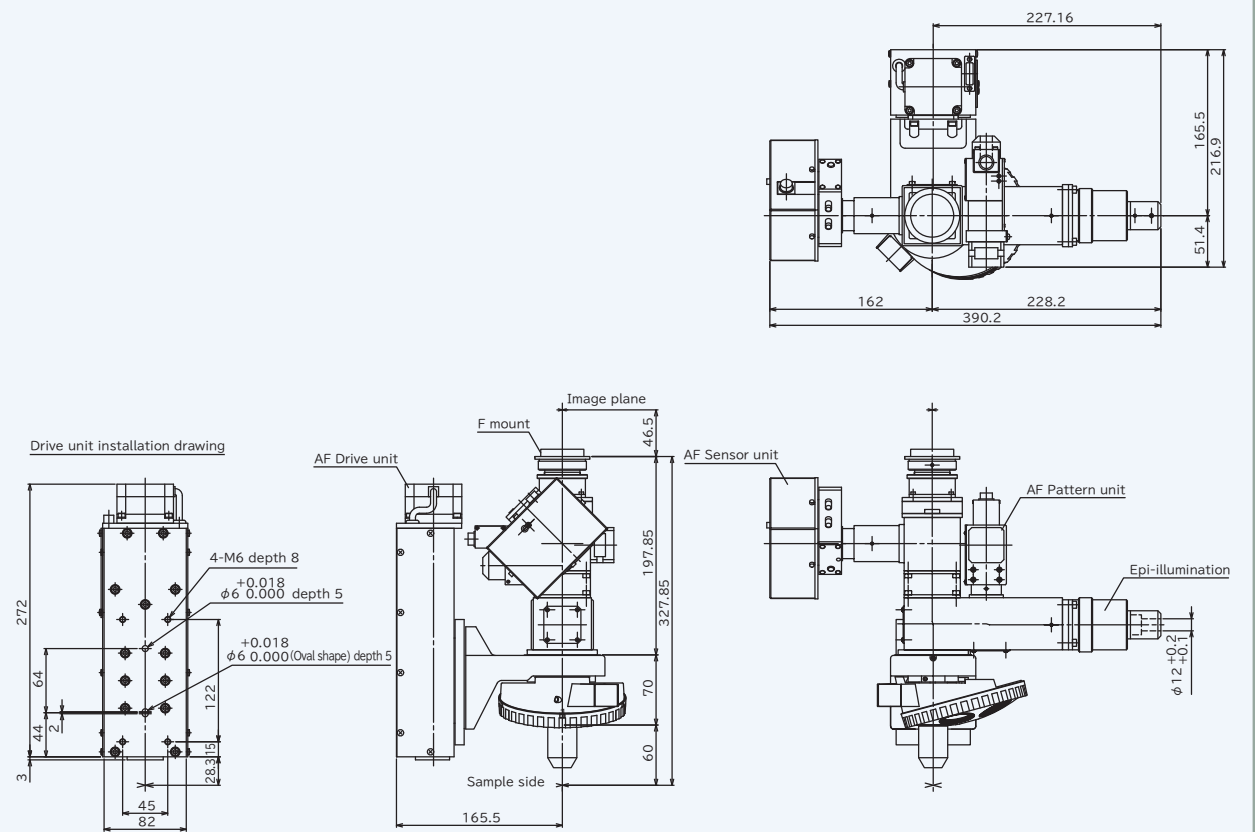
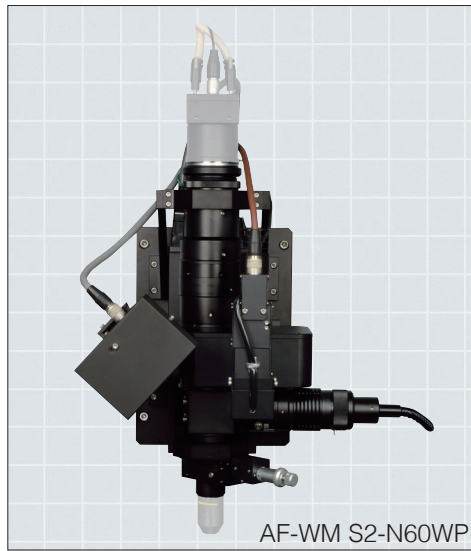
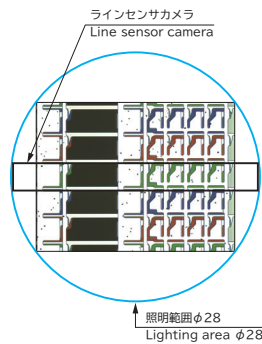
- ① AF-WM is designed to comply with up to φ28 mm or 1.1 inches sensor size, such as C mount area camera and F mount area camera.
- ② This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
- ③ Enables to autofocus on transparent surfaces such as glass and plastics.
- ④ Applicable also to surfaces containing potholes or holes.
- ⑤ Supports a broad range of magnifications from macro to micro.
- ⑥ C mount camera or F mount camera can be used with the dedicated camera mount.
- ⑦ A high-rigidity microscope that is suitable for building into an online machining/measuring apparatus.

*The imaging area vary depending on objective lens and the sample conditions.

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Objective lenses	Nikon, Olympus, Mitutoyo
Camera	C mount camera (Device size 1 inch or less), F mount camera (φ28 mm, equivalent to 1.1 inches)
Drive unit	Model AF-461ZA
	Limiting resolution 0.078 μm
	Travel Standard 4 mm (changeable up to 20 mm)
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)
Focusing	Line sensor method
Focusing conditions (reference value)	Search range ±250 μm (objective lens 20×)
	Search time 0.3 sec
AF Accuracy	Depth of Focus x 1/4

※Maximum cable length 15 m.
 ※Inquire for custom specifications.
 ※For custom specifications, provide us with your objective lenses and camera.
 ※The autofocus accuracy and search time vary depending on the sample conditions, etc.

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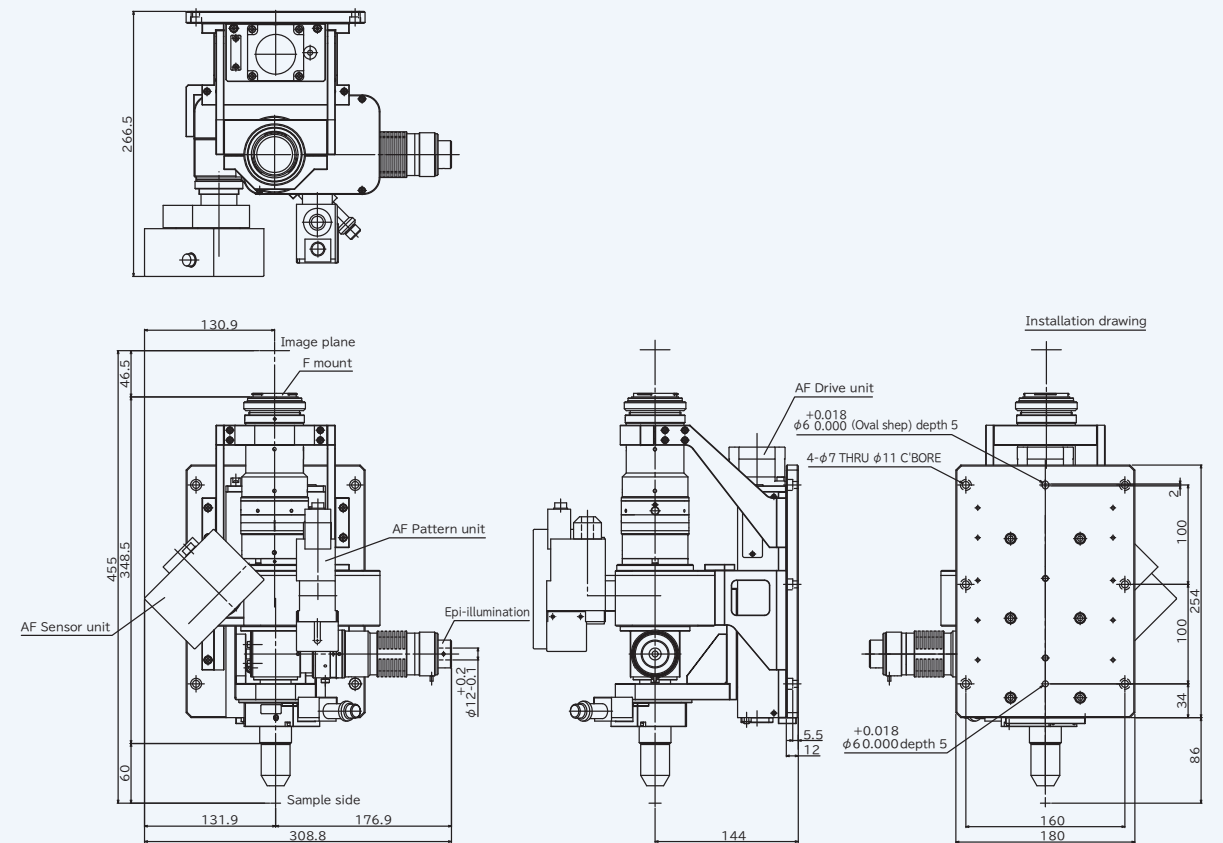
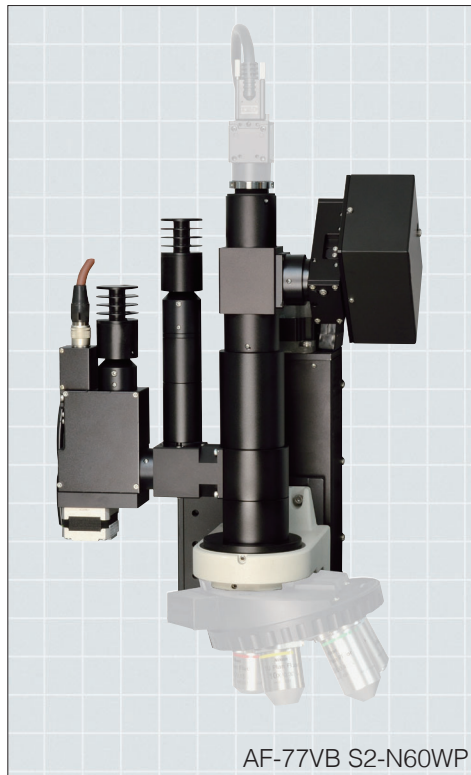
- ① Use of a revolver (or motorized revolver) enables autofocus with objective lens with 2 or more ports.
- ② A compact microscope that can be built into an online machining/measuring apparatus.
- ③ This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
- ④ Enables to autofocus on transparent surfaces such as glass and plastics.
- ⑤ Applicable also to surfaces containing potholes or holes.

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
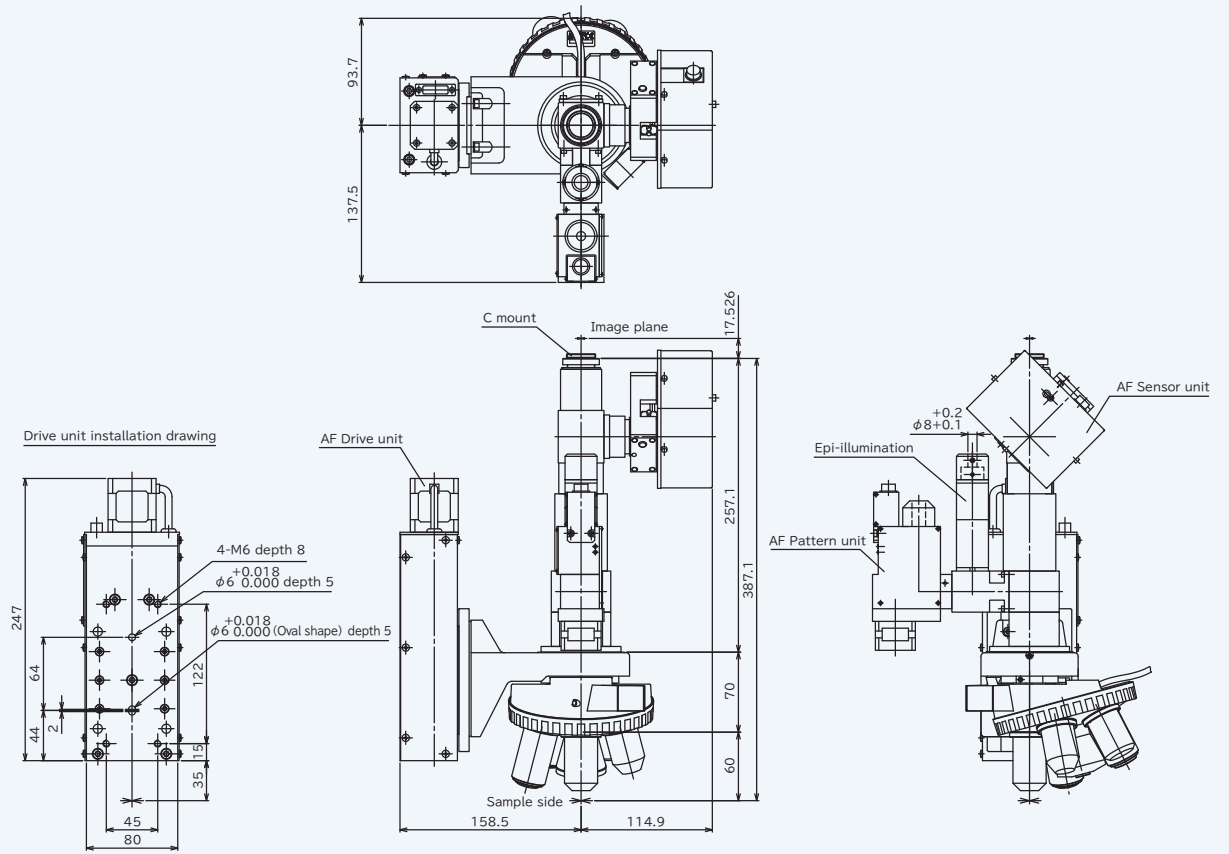
Objective lenses	Nikon, Olympus, Mitutoyo	
Camera	C mount camera (Device size 2/3 inch or less)	
Drive unit	Model	AF-61ZA
	Limiting resolution	0.078 μm
	Travel	Standard 4 mm [changeable up to 10 mm]
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)	
Focusing	Line sensor method	
Focusing conditions (reference value)	Search range	±250 μm (objective lens 20×)
	Search time	0.3 sec
AF Precision	Depth of Focus x 1/4	

※ Maximum cable length 15 m.
 ※ Inquire for custom specifications.
 ※ For custom specifications, provide us with your objective lenses and camera.
 ※ The autofocus accuracy and search time vary depending on the sample conditions, etc.

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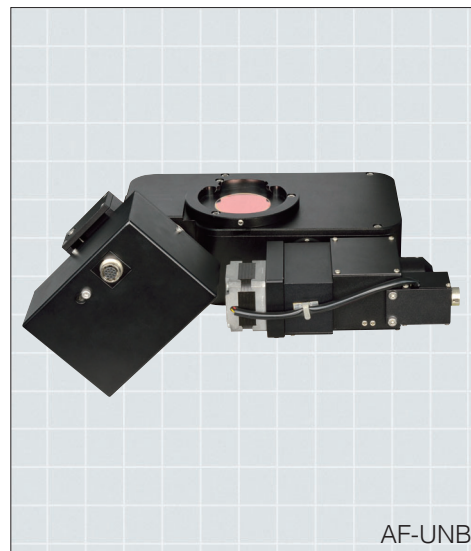- ① Use of an illumination tube and a revolver (or motorized revolver) enables autofocus with objective lens with 2 or more ports.
- ② A compact microscope that can be built into an online machining/measuring apparatus.
- ③ This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
- ④ Enables to autofocus on transparent surfaces such as glass and plastics.
- ⑤ Applicable also to surfaces containing dents or holes.
- ⑥ Easy to install on a commercial microscope and can be used as an autofocus microscope.
- ⑦ Observation methods, BF (Brightfield), DF (Darkfield), Polarizing, and DIC (Differential Interference Contrast) can be chosen depending on the sample.

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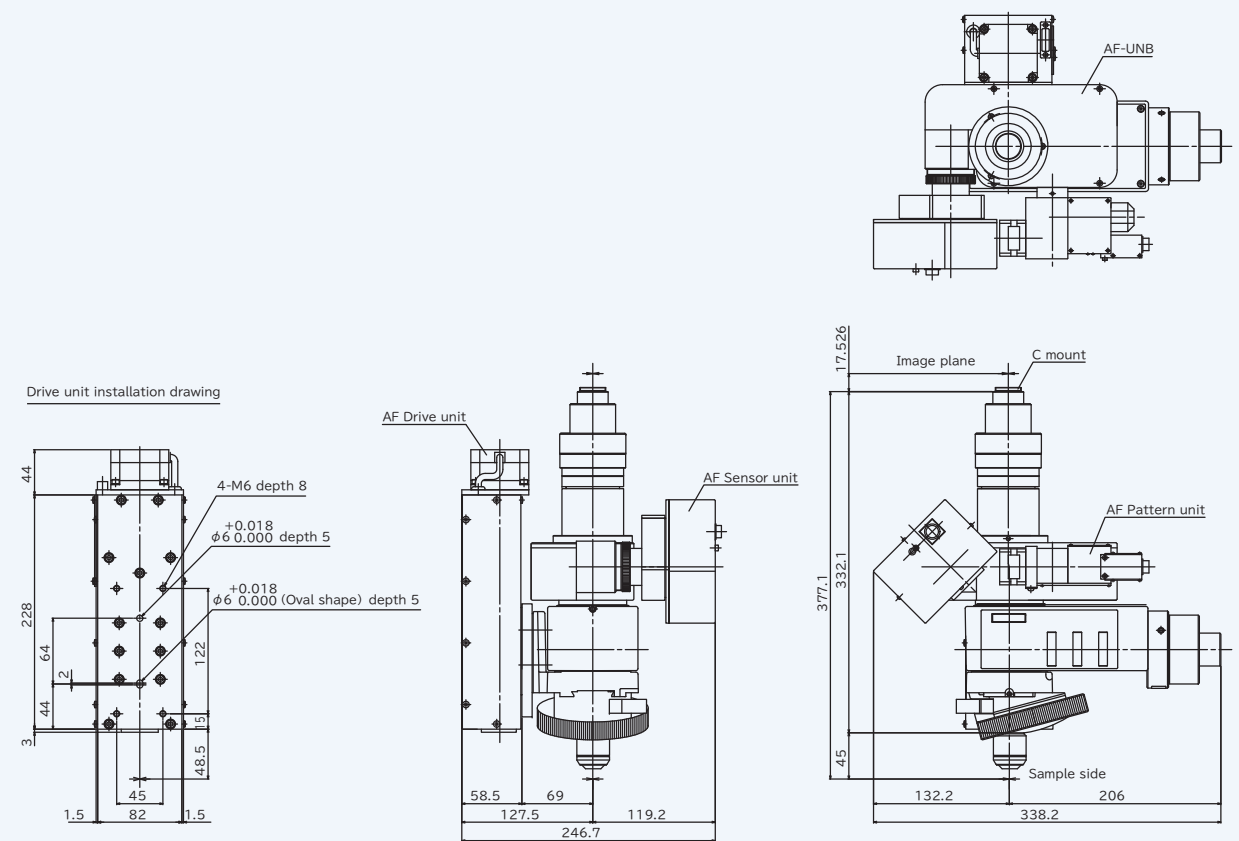
- ① A compact microscope that can be built into an online machining/measuring apparatus.
- ② Use of an illumination tube and a revolver (or motorized revolver) enables autofocus with objective lens with 2 or more ports.
- ③ This microscope is designed to comply with Nikon Corporation CF/IC Series, and CFI60 Series, Olympus Corporation UIS2 Series, and Mitutoyo Corporation objective lenses.
- ④ Enables to autofocus on transparent surfaces such as glass and plastics.
- ⑤ Applicable also to surfaces containing dents or holes.
- ⑥ Easy to install on a commercial microscope and can be used as an autofocus microscope.
- ⑦ Observation methods, BF (Brightfield), DF (Darkfield), Polarizing, and DIC (Differential Interference Contrast) can be chosen depending on the sample.

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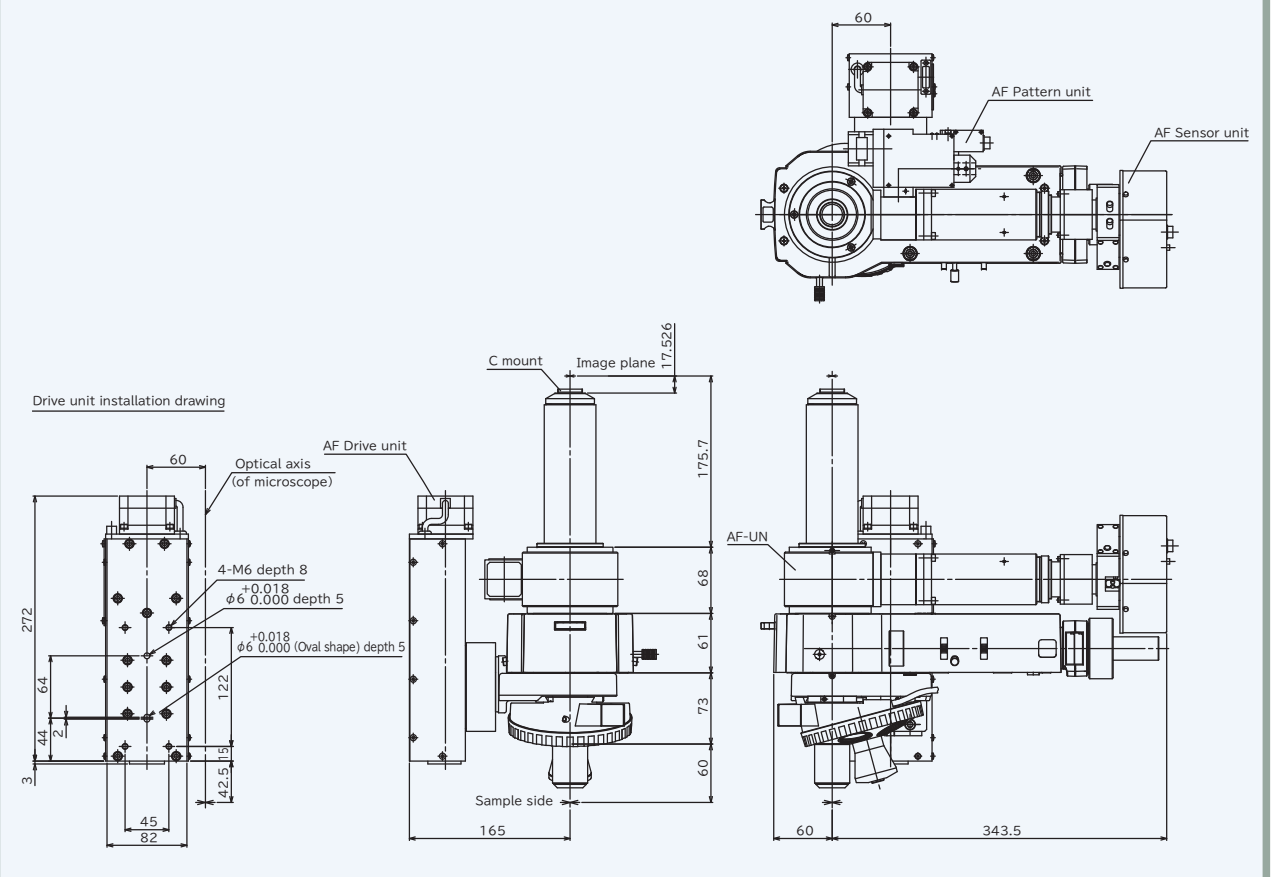
Objective lenses	Nikon, Olympus, Mitutoyo
Camera	C mount camera (Device size 2/3 inch or less)
Drive unit	Model AF-61ZA
	Limiting resolution 0.078 μm
	Travel Standard 4 mm (changeable up to 10 mm)
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)
Focusing	Line sensor method
Focusing conditions (reference value)	Search range ±250 μm (objective lens 20×)
	Search time 0.3 sec
AF Accuracy	Depth of Focus x 1/4

※Maximum cable length 15 m.
 ※Inquire for custom specifications.
 ※For custom specifications, provide us with your objective lenses and camera.
 ※The autofocus accuracy and search time vary depending on the sample conditions, etc.

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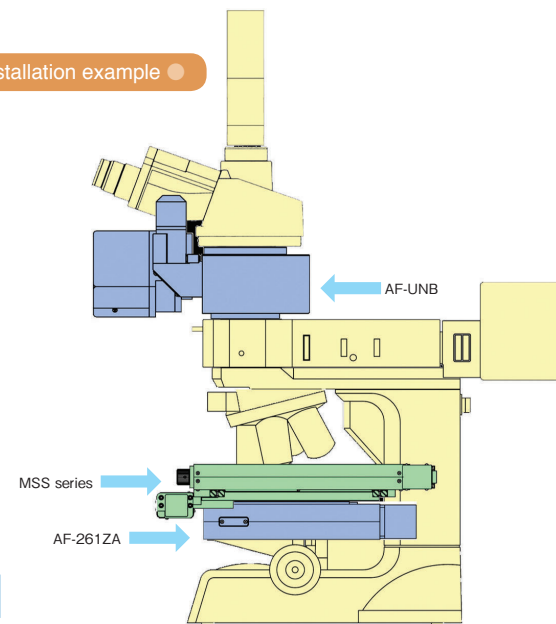
Microscopes	Nikon, Olympus	
Drive unit	Model	AF-261ZA
	Limiting resolution	0.039 μ m
	Travel	2 mm
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)	
Focusing	Line sensor method	
Focusing conditions (reference value)	Search range	\pm 250 μ m (objective lens 20 \times)
	Search time	0.3 sec
AF Precision	Depth of Focus x 1/4	

※ Maximum cable length 15 m.
 ※ Inquire for custom specifications.
 ※ For custom specifications, provide us with your microscope and camera.
 ※ The autofocus accuracy and search time vary depending on the sample conditions, etc.

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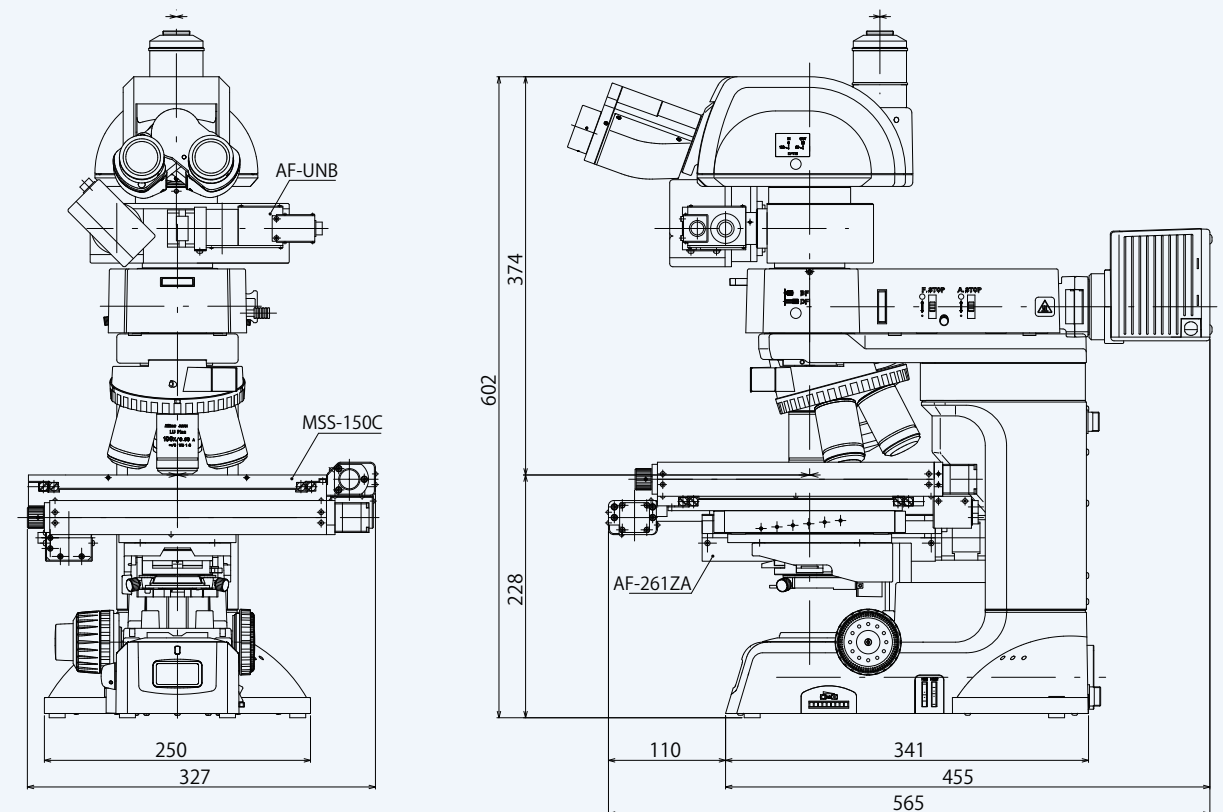
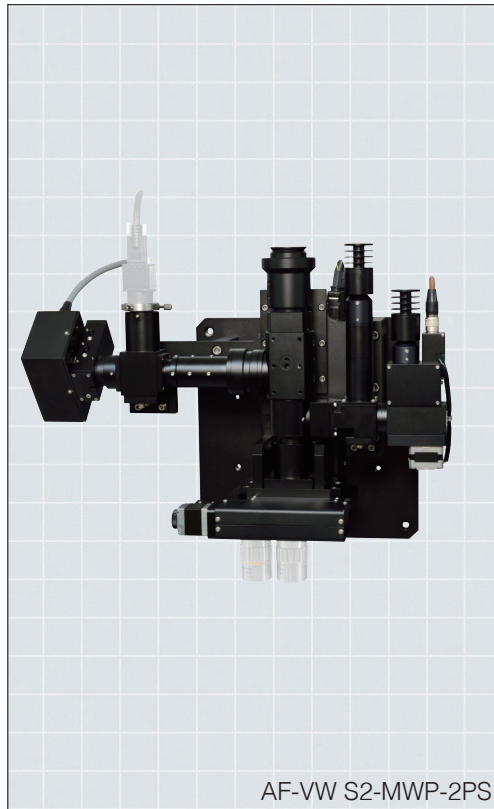
- ① AF-VW is a laser repair system microscope designed for Triple wavelengths (355/532/1064 nm).
- ② Can use more than 2 objective lenses with objective slider. Used for such as microfabrication by laser or camera observations.
- ③ Applicable also to surfaces containing dents or holes.
- ④ Enables to autofocus on transparent surfaces such as glass and plastics.

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Objective lens	Mitutoyo, Shibuya Optical	
Camera	C mount camera (Device size 2/3 inch or less)	
Drive unit	Model	AF-61ZA
	Limiting resolution	0.078 μm
	Travel	Standard 4 mm (changeable up to 10 mm)
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)	
Focusing	Line sensor method	
Focusing conditions (reference value)	Search range	±250 μm (objective lens 20x)
	Search time	0.3 sec
AF Precision	Depth of Focus x 1/4	
Objective Slider	Selectable 2 ports type, 4 ports type or 6 ports type	

* Maximum cable length 15 m.
 * Inquire for custom specifications.
 * For custom specifications, provide us with your microscope and camera.
 * The autofocus accuracy and search time vary depending on the sample conditions, etc.
 * See page 25 for details on AF-OLS Objective Slider.

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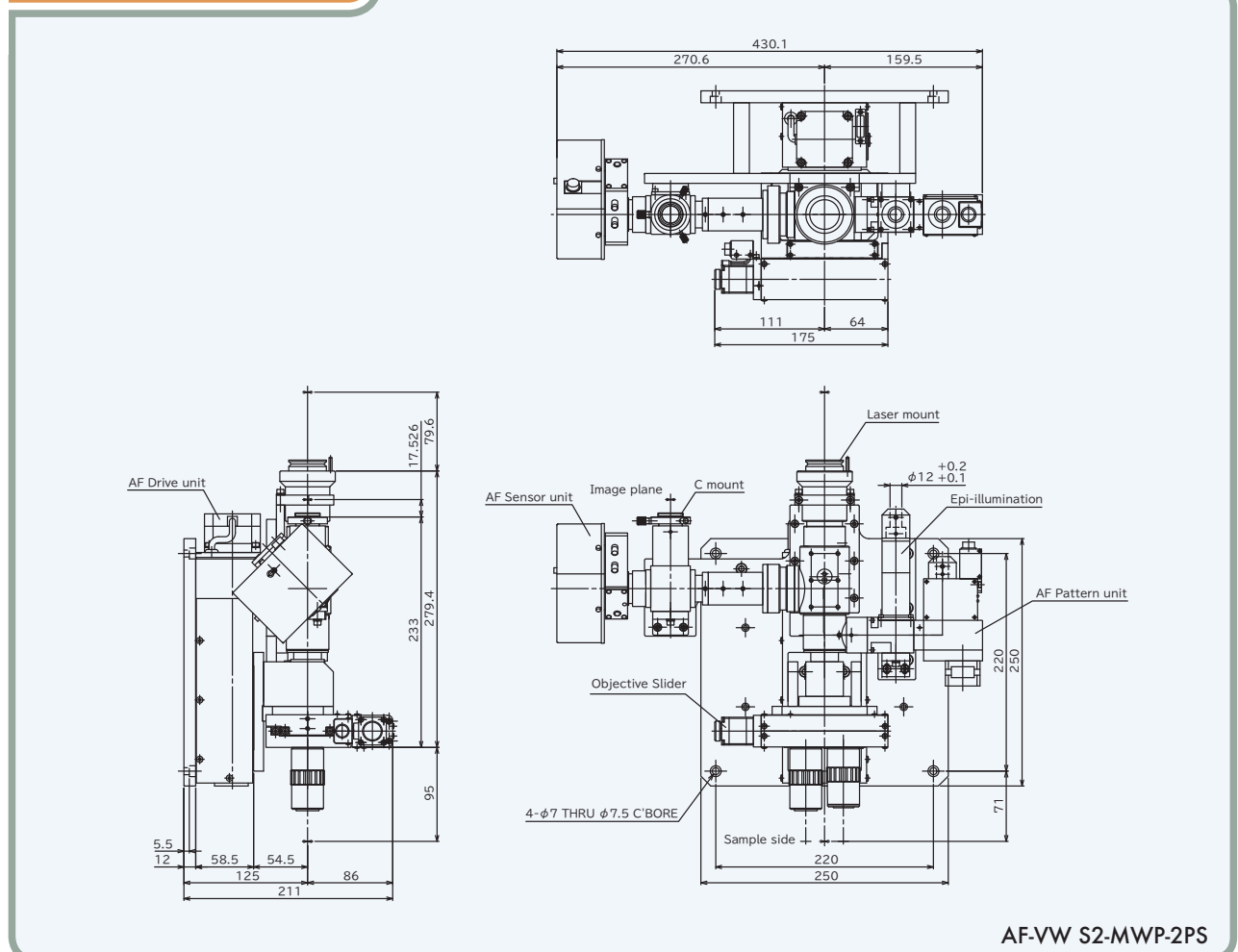
- ① AF-BD is designed to observe simultaneously with Brightfield and Darkfield. No need to change illumination as when using an illumination tube. Brightfield/Darkfield illumination (simultaneous).
- ② Compact microscope with built-in ring illuminator for Darkfield illumination. No need to use an illumination tube.
- ③ Enables to autofocus on transparent surfaces such as glass and plastics.
- ④ Applicable also to surfaces containing potholes or holes.
- ⑤ Capable up to φ28 mm or 1.1 inches image size as when using AF-WF tube system (Wide field type).

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Objective lens	Nikon objective lenses for Brightfield/Darkfield (BD series)	
Camera	C mount camera (Device size 2/3 inch or less)	
Drive unit	Model	AF-61ZA
	Limiting resolution	0.078 μm
	Travel	Standard 4 mm (changeable up to 10 mm)
Cable	Robot cable for sensor, Robot cable for pattern, Robot cable for motor 3 m each (standard)	
Focusing	Line sensor method	
Focusing conditions (reference value)	Search range	±250 μm (objective lens 20×)
	Search time	0.3 sec
AF Precision	Depth of Focus x 1/4	
Ring illuminator for Darkfield	Selectable light guide length 1 m, 2 m, 3 m, or 4 m	

*Maximum cable length 15 m.
 *Inquire for custom specifications.
 *For custom specifications, provide us with your microscope and camera.
 *The autofocus accuracy and search time vary depending on the sample conditions, etc.

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
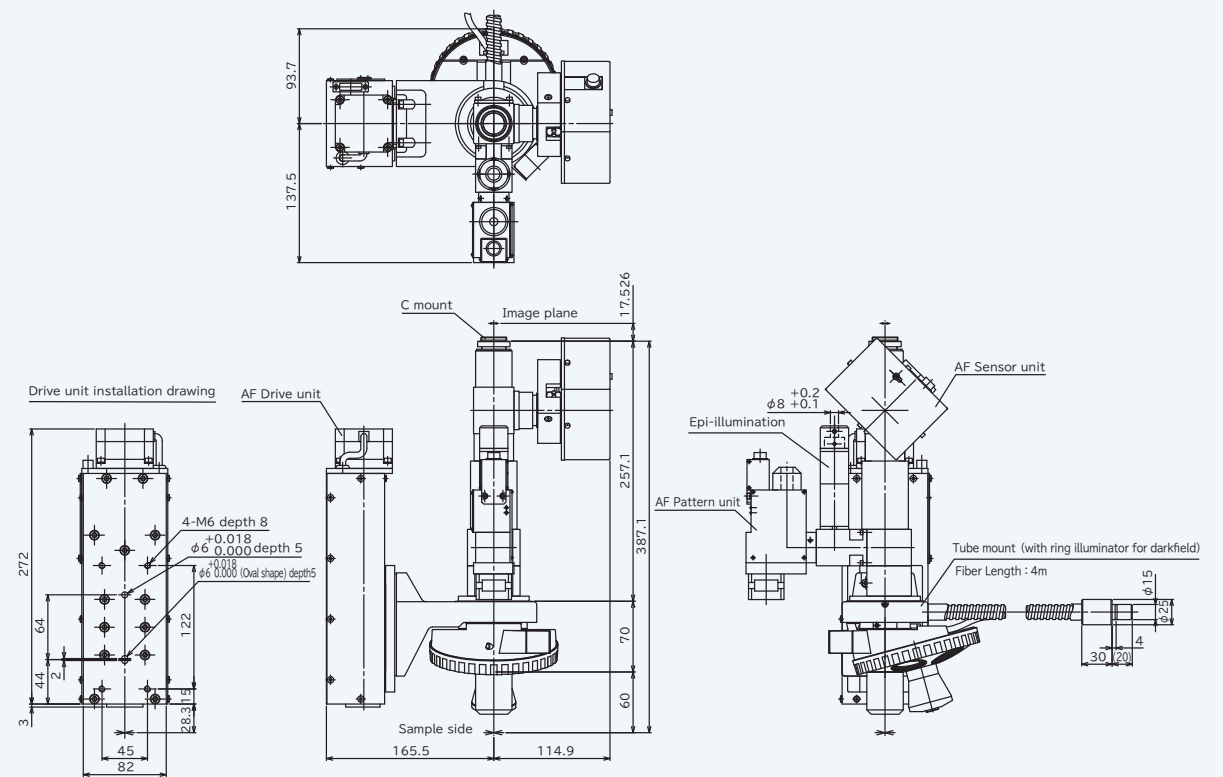
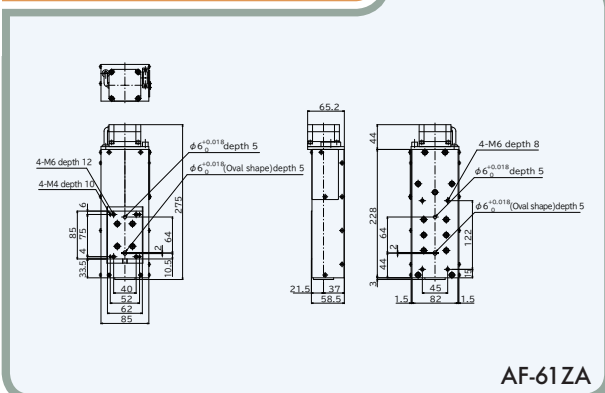
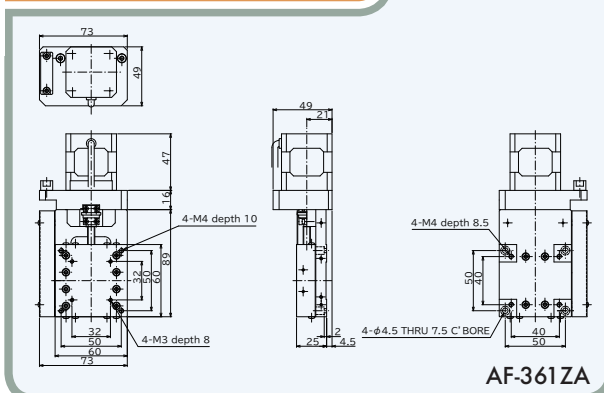
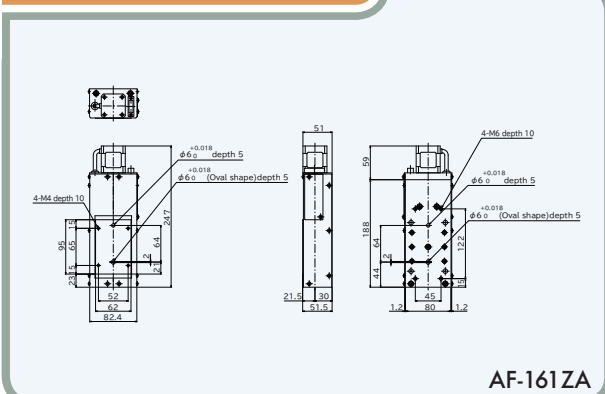
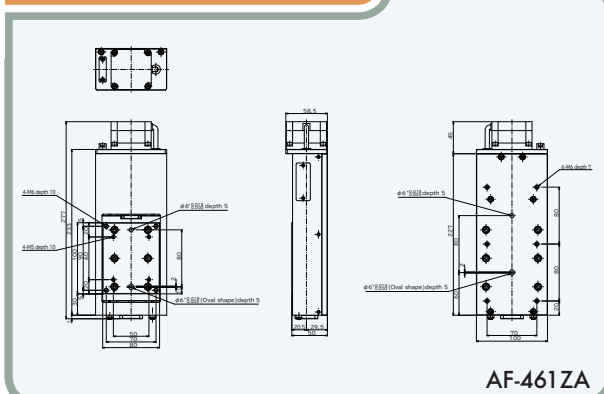
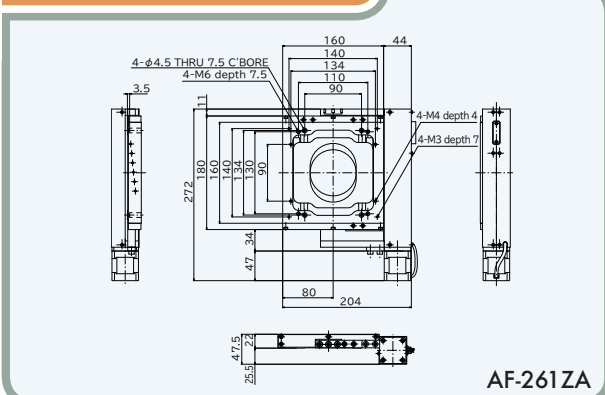
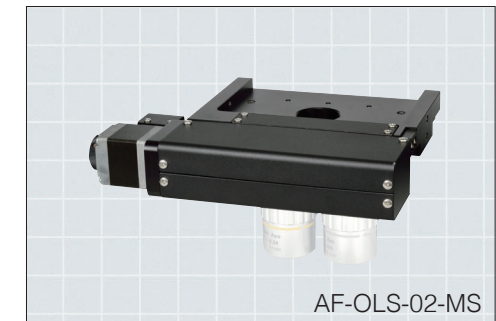
One direction objective slider which enables to switch the objective lens to proper broad range.


- ① AF-OLS, Objective slider has 3 types - 2 ports type (2 objective lenses), 4 ports type (4 objective lenses) and 6 ports type (6 objective lenses).
- ② Can achieve high positioning repeatability with sliding type objective lenses.
- ③ Can be built into a laser repair system and an automatic inspection system, etc.

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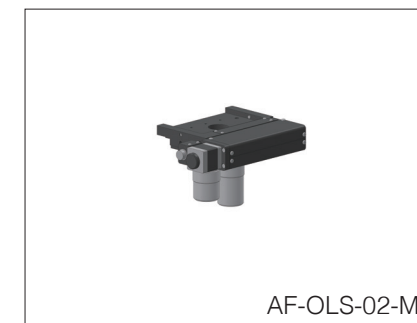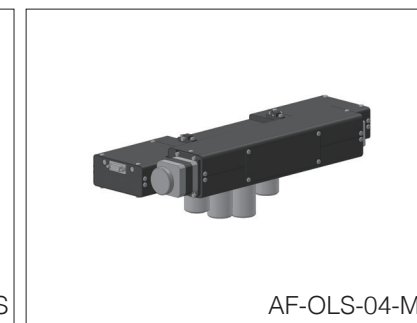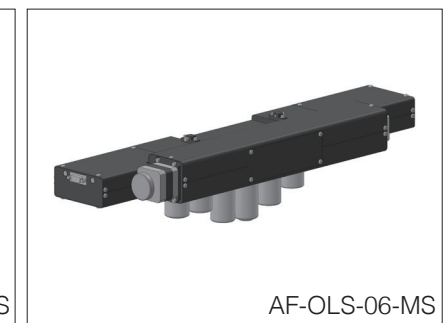
Model	QT-ADL1(K)-AC
Input power	AC 90 to 240 V, 50/60 Hz
Dimensions	W 165 mm x D 140 mm x H 60 mm (not including projections)
Weight	1.6 kg
Control axis	1 axis (with motor driver)
Control	Remote operation (RS-232C)
Communication function	RS-232C

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
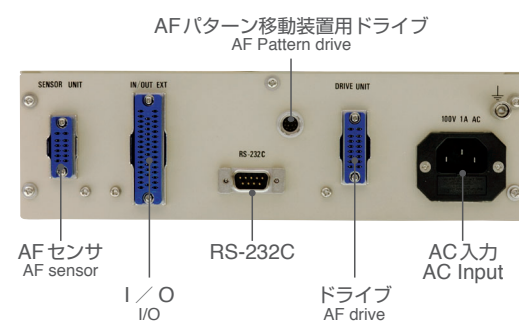
- 1 Controller exclusively for Autofocus.
- 2 Manual operation and remote operation with a PC are possible.
- 3 Special software allows you to configure various settings and controls.

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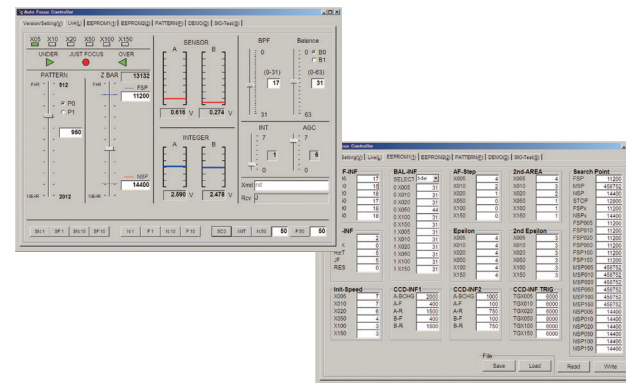
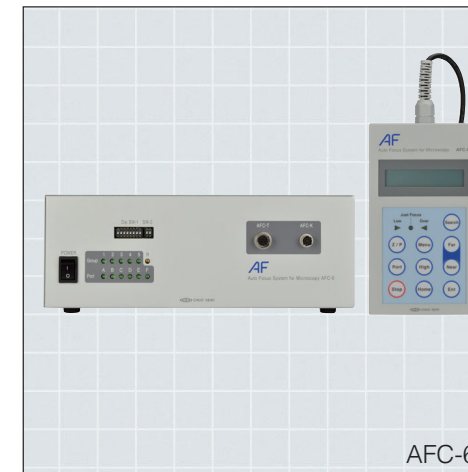
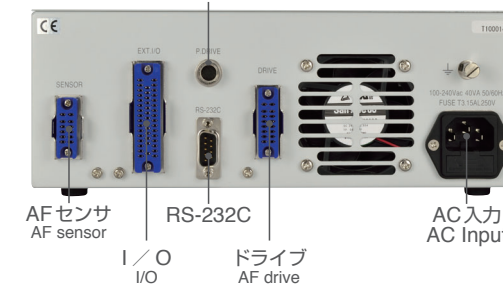
- 1 Controller exclusively for Autofocus.
- 2 Manual operation by Operation box (option) and remote operation with PC are possible.
- 3 Special software allows you to configure various settings and controls.

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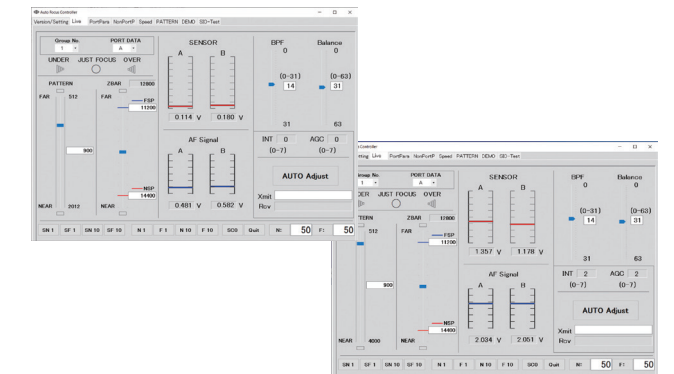General specifications	Power	AC 100 to 240 V, 50/60 Hz
	Current consumption	1A
	Ambient temperature/Humidity	10 to 40°C / 20 to 80% (noncondensing)
	Dimensions	W 254 mm × D 332 mm × H 75 mm (not including projections)
	Weight	3.3 kg
	Option	External operating box (I/O port)
Performance specifications	Target motor	2 phase stepper motor, 1 A/phase or less (depending on the conditions)
	Control	Stepped shift, return to origin, port operations, remote operations (option)
	Drive frequency	Max. 20 kHz
	Return to origin	Machine zero, offset zero
	Over run	Limit switch
	Communication function	RS-232C, I/O

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General specifications	Power	AC 100 to 240 V, 50/60 Hz
	Current consumption	1A
	Ambient temperature/Humidity	10 to 40°C / 20 to 80% (noncondensing)
	Dimensions	W 250 mm × D 280 mm × H 90 mm (not including projections)
	Weight	3.3 kg
	Option	Operation box (AFC-K)
Performance specifications	Target motor	2 phase stepper motor, 1 A/phase or less (depending on the conditions)
	Control	Stepped shift, return to origin, port operations, remote operations (option)
	Drive frequency	Max. 20 kHz
	Return to origin	Machine zero, offset zero
	Over run	Limit switch
	Communication function	RS-232C, I/O